


Focus&Zoom Operation Guide for Item 2319418/ BW1008-500X



1. Fix a viewing distance(0-25mm) between microscope and object, try to make microscope much closer to object.



2. turn the focus knob to the far left(small magnification side), make it to be minimum; then turn the knob to the right(big magnification side) slowly and slightly, until appeared clean image under a smaller magnification, and then go on turning, until appeared another clean image under a bigger magnification.

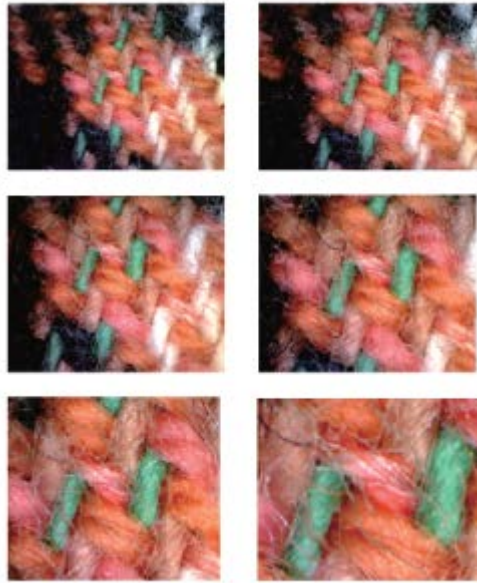

3. the final step is to press the “zoom” button, to do zooming: when you press “zoom” button one time, the object will magnified 1.3 times, press twice, it will magnified another 1.3 times, go on pressing until magnified the maximum, at this time if go on pressing, the image will resume the original size, you can process it circularly.