



Introduction

This is a smart-handly size Digital Microscope! It can easily link to your PC and applicated well in different fieds :
 A Electronic Filed: PCB and components inspection
 B Material: Powder, Material surface inspection
 C Beauty shop: Hair & Skin inspection
 D Education: Find a small world
 Based on PC supports, this Digital-Microscope can do many powerful works: It can provide more friendly functions: Snap a digital photo, Recording Video and its compact size makes your observation easily. Digi-Microscope will support users do their works more easily and never miss a detail.

Parameters And Features


- * 300K 480K 1300K Pixels Resolution
- * High performance CMOS sensors
- * Window capture size: 640X480
- * Video mode: 24bit true color
- * Interface: USB 2.0 / USB 1.1
- * Transmission rate: 640X480 30 f/s
- * Focus Range: 5mm-45mm(FFL)
- * Buld-in image compression function
- * Automatic brightness adjustment
- * Automatic color compensation
- * Manual focus

System requirements


- * IBM or compatible PC with USB port
- * Pentium 200 or higher CPU
- * Windows 2000 / XP / Vista / Win 7
- * Over 200MB free harddisc space
- * CD-ROM drive
- * 32MB or higher memory
- * Support DirectX VGA card

Driver installation

1. Insert Driver/Application disc into CD-ROM drive and open driver directory. Double click Driver-setup\Setup.exe and the system will automatically run necessary files and display a window as following

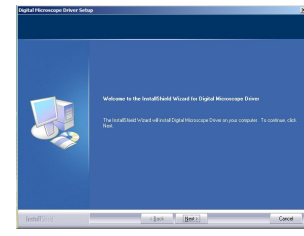

Click "Next"

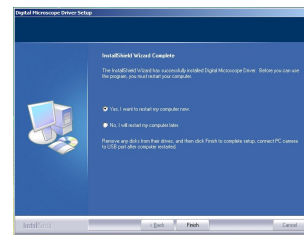


Click "Finish"


2. Plug-into USB port in PC, the system will find it and automatically complete the installation



When installation is completed, you will find Amcap from Programs

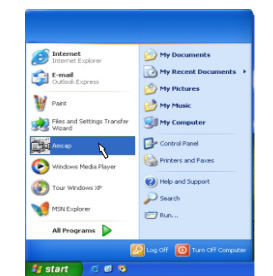

3. Run Amcap.exe and you will get the picture on your PC display.

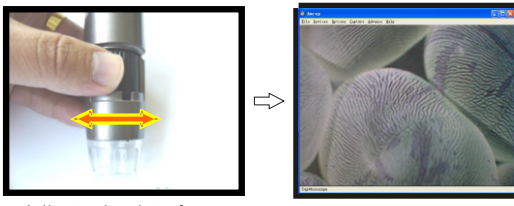


Adjust wheel to focus a clear image

4. Press snap button to make a quick start amcap provide many useful program for photo treatment.




How To Make A Perfect Photo

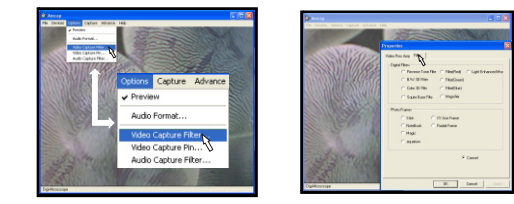


A: Select Video Capture Filter B: Click "Filter"

Digital Filters

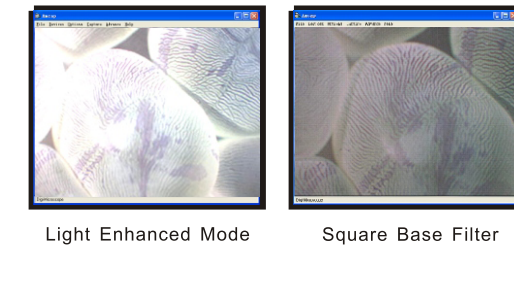

Light Enhanced Mode Square Base Filter

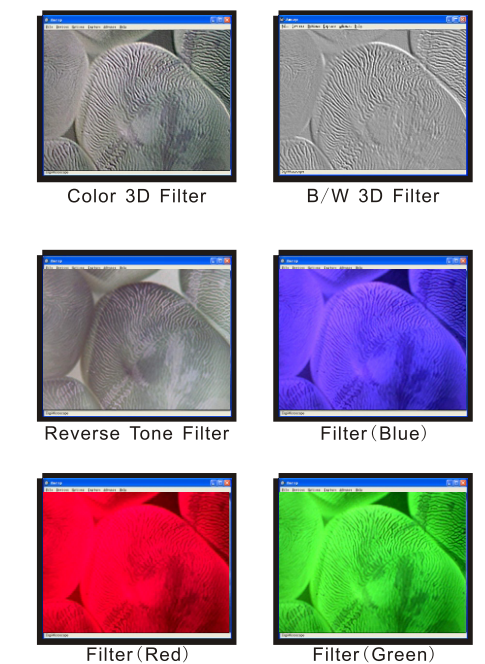

Color 3D Filter B/W 3D Filter
 Reverse Tone Filter Filter (Blue)
 Filter (Red) Filter (Green)

Photo Frame



Photo Frame 1 Photo Frame 2


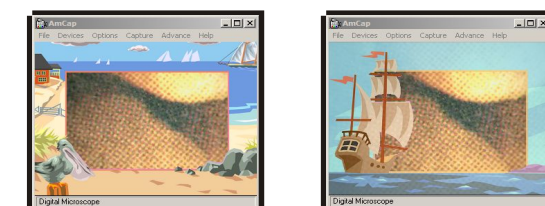


Photo Frame 3 Photo Frame 4

Select Image Resolution (Size)

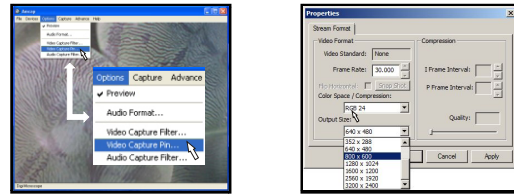

A: Select Video Capture Pin B: Choose the size you need

Caution

- * Avoid contact to the lens with hand, sharp or coarse stuff; use clean and soft cloth for cleaning
- * Do not use it in hot, cold, dusty or humid circumstances
- * Avoid falling as much as possible. Severe falling may cause damages to the lens, malfunction or scratches to the surface

Warranty

- * This product is under warranty of non-human factors:
 - A. Free replacement within three months from purchase date
 - B. Free charge for repair within one year from purchase date
- * For any problems concerning our products you purchase, please send them back to local sales agent with effective note